REMOTE CALIBRATION UNIT





USER MANUAL

۲

Th/Ar Ref.On

Flat Ref.Off

.

Remote

RCL

February 2015 – English Revision v0.3

Baader Planetarium GmbH

Zur Sternwarte D - 82291 Mammendorf Tel.: +49 (0) 8145 - 80 89-0 Fax: +49 (0) 8145 - 80 89-105 www.baader-planetarium.de

Remote **C**alibration **U**nit Professional Set for BACHES

User Manual

Thank you for purchasing this Baader Planetarium product. To get the most out of your RCU calibration set, please read this instruction manual thoroughly before starting to work with the RCU unit. Keep this user manual available for reference and visit the Baader Planetarium website: http://www.baader-planetarium.de for new information about the product.



Unpacking

As you unpack the RCU, check the list below and make sure that all items shown above are included. If anything is missing, contact your original retailer.

Packing list for the standard model

- Remote Calibration Unit
- Power cord for EU plug socket
- Fiber docking port for BACHES (already mounted on BACHES if RCU was ordered together with BACHES)
- Glass fiber to connect RCU with spectrograph
- Power cord for the motor in BACHES
- Watertight, shock-resistent ABS flight-case with machined compartments to ensure maximum transport safety (not shown)
- USB stick with software and contend
- Manual



Part description

- 1. Glass fiber coupler
- 2. Power connector for motor
- 3. Three position switch for OFF, Th/Ar ON, or flat-field ON
- 4. Two position switch to enable (Ref.On) or disable (Ref.OFF) the coupling calibration mirror in BACHES
- 5. Two position switch to select remote and manual operation



Part description (continued)

- 6. ON/OFF switch
- 7. 220V Connector
- 8. 2A Fuse
- 9. Remote Control Unit panel

	8
1.1 The main features of the RCU	8
1.2 Handling advice	9
2 QUICK SETUP	10
2.1 Getting started	10
2.2 Installing the fiber coupler to your BACHES	11
2.3 Attaching the fiber	12
3 STARTING THE RCU	13
3.1 Activate the RCU	13
3.2 Activate the Halogen lamp	13
3.3 Activate the Thorium-Argon lamp	14
4 RCU REMOTE CONTROL	15
4.1 Activate remote control	15
4.2 Remote control via the RS232 serial port	15
•	17
4.3 Remote control via LAN and web interface	···· /
4.3 Remote control via LAN and web interface	
4.3 Remote control via LAN and web interface	
 4.3 Remote control via LAN and web interface 4.3.1 Ethernet LAN switch on the RCU 4.3.2 Status LED 4.3.3 Manual IP addresses 	17
 4.3 Remote control via LAN and web interface	17 17 17 18 19
 4.3 Remote control via LAN and web interface	17 17 18 19 19
 4.3 Remote control via LAN and web interface 4.3.1 Ethernet LAN switch on the RCU 4.3.2 Status LED 4.3.3 Manual IP addresses 4.3.4 IP addresses via DHCP 4.3.5 Configuring by Software 4.3.6 Web interface 	
 4.3 Remote control via LAN and web interface 4.3.1 Ethernet LAN switch on the RCU 4.3.2 Status LED 4.3.3 Manual IP addresses 4.3.4 IP addresses via DHCP 4.3.5 Configuring by Software 4.3.6 Web interface 	
 4.3 Remote control via LAN and web interface 4.3.1 Ethernet LAN switch on the RCU 4.3.2 Status LED 4.3.3 Manual IP addresses 4.3.4 IP addresses via DHCP 4.3.5 Configuring by Software 4.3.6 Web interface 5 SPECIFICATIONS 6 USEFUL ACCESSORIES	

1 Introduction

1.1 The main features of the RCU

"RCU" is an acronym for **R**emote **C**alibration **U**nit. The RCU is a multifunctional spectrograph calibration tool mainly designed for use with the BACHES spectrograph. It contains a Thorium-Argon lamp and a filtered Halogen flat-field lamp to work as a standalone calibration unit.

RCU's main features are:

- Thorium-Argon calibration lamp incl. a speciallydesigned power supply
- Lens system to inject light beam from Th/Ar spectral calibration lamp into the glass fiber with the highest efficiency
- Blue filtered Halogen lamp for the purpose of flatfielding
- Switchable 12V power output for the coupling flipmirror on BACHES
- Remote unit with LAN or RS232 connection
- Factory tested and aligned

1.2 Handling advice

The RCU includes a calibrated optical system to feed the light of either the Th/Ar spectral calibration lamp or the Halogen flat-field lamp with the highest efficiency possible into the fiber. Please handle the RCU with care. Hard shocks may destroy the optical alignment and reduce the efficiency dramatically. The Th/Ar lamp is well protected in a fixed unbreakable encasement. However the lamp can break if the device gets a hard blow or hits the ground.

The amount of Thoriumoxide in the bulb is below the saftey threshold limits. Nonetheless, it is not healthy to inhale the gases if the lamp breaks, just as if an energy-saving lamp were broken. We suggest opening all the windows exposing the RCU to fresh air for some hours. The surface of the unit should be wiped with a clean soft cloth or soft paper tissue, subsequently discarded.

Please carefully read the material safety data sheet¹ provided on the USB stick. If the device is misaligned or the lamp is broken, faulty or worn out, the RCU must be returned to Baader Planetarium GmbH for replacing the lamp and realignment.

Transport regulations do not classify this product as dangerous. It is mandatory not to discard the lamp to avoid environmental contamination. Disposing such lamps in local deposits or household garbage is strictly forbidden. Baader Planetarium GmbH does not take any responsibilities in such case. Broken or used lamps must be returned to Baader Planetarium GmbH for disposal.

¹ https://www.edwardsvacuum.com/uploadedFiles/Resource/MSDS/MSDS_German/P120-10-000-882C.pdf

The RCU has a regulated power supply adjusted to the lamp current suggested by the manufacturer. Based on the current, the lamp has a lifetime of 500 hours. It may degrade during that time, leading to a longer integration time even if the RCU is still calibrated. Degradation is normal and must be accepted. In case the user forgets, the RCU has a built-in timer to turn off the Th/Ar lamp after 20 minutes. This feature is included to extend the lamp's lifetime.

2 Quick Setup

2.1 Getting started



Put the RCU on a firm support. Plug in the power cord into a wall socket.

The RCU needs 200V-240V mains voltage. If your national power grid provides 110V, the RCU be run with a 500W step-up transformer.



Place your spectrograph close to the RCU.

2.2 Installing the fiber coupler to your BACHES

This has already been done if the RCU was ordered together with BACHES. If so, skip ahead to section 2.3.



Remove the cover installed on your BACHES. Mount the cable-holding bracket that comes with the RCU.



Your BACHES should now look like this.



Gently fasten the fiber coupler to BACHES head using the supplied screws. The fiber coupler may need some alignment to work properly; how to do that is explained later on in this manual.

2.3 Attaching the fiber



Attach the fiber to the fiber coupler. Take care to not scratch the polished front window of the fiber while plugging it in.

Attach the power cord to the motor cover.

Secure the fiber and the power cord to the holding bracket.



Attach the fiber and the power cord to the **front connectors** of the RCU. Do not try to plug on the RS232 socket at the rear side of the RCU. You are now ready to start working with the RCU.

3 Starting the RCU

3.1 Activate the RCU



Shift the right switch to **Manual** and the middle switch to **Ref. On**. This will flip the coupling mirror in the BACHES head into the calibration position. For accurate calibration, ONLY calibration lamp light may illuminate the BACHES'

slit. When the reference switch is ON, be sure to cover all openings of the BACHES head with their dust caps to exclude any stray light that will interfer with calibration.

3.2 Activate the Halogen lamp



Shift the left switch to **Flat** to enable the flat-field lamp. Attach an eyepiece or camera to BACHES per its instructions manual.

You should see a flat-field BACHES echelle spectrum as shown to the left. Usual integration time may be five to ten seconds.

3.3 Activate the Thorium-Argon lamp



Switch to the **Th/Ar** lamp. The integration time for Th/Ar is usually a little shorter then for the flat-field lamp.

Every Th/Ar lamp needs some time to stabilize. To achieve very

precise calibration, turn on the Th/Ar lamp at least five to ten minutes beforehand. This will consistently give the same intensity level.

While the Th/Ar lamp is switched on, do not look directly at the fiber output. UV radiation from the lamp may harm your eyesight.



Thorium-Argon spectrum obtained with a Nikon D7100 DSLR camera, exposure time 10s at ISO 1000

4 RCU Remote control

4.1 Activate remote control



Shift the right switch to **Remote** to enable remote control

4.2 Remote control via the RS232 serial port



To run the device via RS232 serial port, you need an optional RS232 adapter cable. Please contact Baader Planetarium. You also need a terminal program such as the Windows application "HyperTerminal". Unfortunately, this tool is no longer implemented in Windows 7, and its status in Windows 8 is unknown. You may download HyperTerminal for free from: <u>http://itler.net/windows-7hyperterminal-so-bekommt-man-es-zurueck/</u>

- 1. Download/install a terminal program, e.g. HyperTerminal
- 2. Connect your PC to the RCU via an RS232 cable
- 3. Switch the RCU ON. When the STATUS LED glows green, choose the COM port to which the RCU is attached. To know which COM port that is, check the device manager of your PC. Please note that a test with Windows 7 and HyperTerminal worked very well even if the status LED stayed red.

Terminal program settings

Bits per second	115200
Databits	8
Parity	None
Stoppbits	1
Flowcontrol	None

If you do not use HyperTerminal, make sure that your terminal program supports VT100 commands.

🇞 RCU-RS232 - HyperTerminal		
File Edit View Call Transfer Help		
Expert Power Control 1002 v2.0.1		
<pre><1> OFF - ThAr-Lampe <2> OFF - Halogenlampe <3> OFF - Einspiegeln</pre>		
<pre><4> UFF - Power Port 4</pre>		
press <1><4> to toggle port 14 press <z> for configuration menu press <esc> to logout</esc></z>		
-		

If the connection is successful, the RCU replies with a login screen. Now the power ports can be switched ON and OFF by typed commands from the PC. An explanation of available commands will be shown as well.

4.3 Remote control via LAN and web interface

4.3.1 Ethernet LAN switch on the RCU

Connect your RCU to your PC by an Ethernet LAN patchwork cable via a 100/1000 Mbit switch. Direct connection might not work unless you use a LAN crossover cable (cables are not supplied).



4.3.2 Status LED

Red	Not connected to Ethernet
Orange	Connected to Ethernet, waiting for TCP/IP settings
Green	Connected to Ethernet, TCP/IP settings finished
Blinking	Unit in Bootloader modus

4.3.3 Manual IP addresses

IP address	192.168.0.2
Subnet	255.255.255.0
To enter the RCU Web interface simply change	your PC Network
settings to:	

IP address	192.168.0.1
Subnet	255.255.255.0

Please follow the instruction in the attachment, if you do not know how to change IP Settings of your PC.

Later on, you can change the IP Settings of the RCU directly in the web interface.

Expert Power Control 1002 - C +	
(I192.168.0.2/config_ip.html	マ 🕑 - Google 👂 🏠 自 🦊 🏫 🛐 🚍
Control Panel Configuration Logout Power Ports - IP Address	· IP ACL · HTTP · SNMP · Syslog
Configurat	ion - IP Address
Hostname:	RCU
• IP Address: • Netmask:	255.255.255.0
• Gateway:	192.168.0.2
• Use DHCP	● yes [®] no
	Арріу
Expert Pov	ver Control 1002 - v.2.0.1

4.3.4 IP addresses via DHCP

If you integrate the RCU into an existing Ethernet with DHCP Server, the RCU the DHCP server will allocate an idle IP address. Please check your DHCP Server for the IP address allocated.

You may want to allocate a unique MAC address in your DHCP Server. The MAC address of your individual RCU unit is found on its serial number documentation page.

4.3.5 Configuring by Software

If your RCU can not be addressed either initially or after changing IP settings in the web interface, please use the supplied software **GBL_Conf.exe**, found on the provided USB stick.

Start the program and start search > All Devices . This will show your RCU. You can now change the settings according to your needs.

4.3.6 Web interface

Start any browser and enter the RCU's IP address 192.168.0.2 or any other IP address you allocated to the RCU during the above start-up procedure.

Now you can fully control your RCU.

Expert Power Control 1002 - R +	-				_ 0	x
	⊽ C ⁱ	<mark>8</mark> ₹ Google	오 🚖 🗎	÷	f 3	≡
Control Panel Configuration Logout						
	Switching · Batc	hmode				
Th/Ar L	amp	On	Off Reset			
📴 Flat Lar	np	On	Off Reset			
FF Referer	nce mo	de On	Off Reset			
💴No de	vice	On	Off Reset			
	(autorefresh: 1	0s)				
Expe	rt Power Control 1	002 - v.2.0.1				

You can switch the RCU ON and OFF simply by typing the buttons of the LOGIN window. Then you can simply switch the devices ON and OFF as shown above.

Be aware that Th/Ar lamp and flat-field lamp can both be switched ON at the same time when using the web interface. This will not damage the RCU, but the image showing both spectra is useless.

Do not forget to switch the RCU to "reference mode On" before starting the calibration.

If you want to use other settings like HTTP access, SNMP or Syslog please refer to the detailed instructions of the remote control unit (PDF on USB stick). We supply a copy of the original manual with each RCU.

5 Specifications

Efficiency ratios for both lamps vary between 1:1 and 1:1.5 depending on the individual unit. We cannot guarantee for a specific number.

Dimensions	345 x 217 x 128 mm
Weight	5,8 kg
Power	Input 200 – 240V / AC, 2A
	50 / 60 Hz

6 Useful accessories



The rack holder mounting plates (#2458623), including cage nuts, are for mounting the RCU either forwards or backwards in a 19" server rack.



The telescope mounting bars (#2458622) are made of stainless steel, and provide a more stable mounting platform than the usual rubber bumpers if the RCU is used outside the laboratory.

7 Troubleshooting

Problem	Suggestion
The Th/Ar lamp suddenly quit working.	The automatic 20 minute time-out may have turned it off. Try turning the lamp OFF and back ON again.
I don't get any light into the BACHES, either Th/Ar or flat- field.	Your optical fiber may not be aligned properly. Please recheck the alignment. It is also possible that the fiber is damaged. Before reinstalling it, see if light is visible at the polished face of the fiber when one of the lamps is on; if not, the fiber must be replaced. Attention: please be aware that part of the light from the Th/Ar lamp consists harmful UV radiation. If possible use only the flat field lamp for visual checks or use sunglasses with UV-blocker for checking Th/Ar light. Do not stare unlimited into the beam.
My flat-field and calibration spectra are very dim.	This may also be a fiber alignment problem. Please recheck the alignment.
The flat-field light is fine, but my Th/Ar is very dim now.	How many hours do you have on the lamp? We suggest keeping a lamp time log. If it is approaching 500 hours, a new lamp may be in order.
My calibrations are not accurate. Everything looks normal, but they don't work.	You may have some light leakage. Are all the dust caps on the BACHES head in place and in good con- dition? If any have cracked or no longer fit snugly, they should be replaced. If you are using the web interface, make certain that only the Th/Ar lamp is ON, and the flat-field lamp is OFF. If you are using a camera to capture the spectrum, make sure that no light leakage is coming in via the camera viewfinder (cover it).
My unit is not working at all. It has power to it.	Are you in a country with 110V power? If so, do you have the appropriate step-up transformer between the wall socket and the RCU?
My Th/Ar spectrum has changed on me a lot, and it no longer looks like what I remember.	The bulb may be compromised by gas leaking in or out of it. Was the unit accidentally bumped hard just before this started? If so, the lamp may need to be replaced.

COPYRIGHT NOTICE

This document is proprietary and for use of the intended recipient only. Baader Planetarium GmbH reserves all rights. All information contained in this document is subject to Baader Planetarium's copyright.

Except for brief quotation in critical articles or reviews, no reproduction of this manual, in any form, in whole or in part, may be made without written authorization from Baader Planetarium GmbH.

Any unauthorized copying, any publishing of its content in the internet or intranet, any use of this information by third parties, and/or its dissemination to third parties, without the expressed written consent of Baader Planetarium GmbH, is a serious infringement.

Under international copyright laws, any reproduction or dissemination of this document by the recipient is expressly prohibited. Any transfer of this document, in original or copied form, to competitors or other third parties violates applicable public copyright laws. Violations of copyright or public procurement laws will be prosecuted under the law –

© Baader Planetarium GmbH, 2015

REMOTE CALIBRATION UNIT

RCU-Unit with BACHES Spectrograph

www.baader-planetarium.de

DER PLANE



BAA

Zur Sternwarte • D-82291 Mammendorf • Tel. +49 (0) 8145 / 8089-0 • Fax +49 (0) 8145 / 8089-105 Baader-Planetarium.de • kontakt@baader-planetarium.de • Celestron-Deutschland.de

 Δ

MB